

Thermostatic device “Set up” mode

1. Switch off the power supply .
2. Press and hold the buttons SET and Start simultaneously and hold them pressed.
3. Switch on the power supply.
4. In 2 seconds inscription SET appears on the display, accompanied by double sound signal.
5. Release the buttons. Now the display shows 40.0 - which is the initially set temperature. By pressing the button SET you can increase the temperature to 65 C. Next pressing returns it to 40 degrees. Press the button START to save the set temperature.
6. Setting up “Cycle 1” time (Indicator Timer 1 lights) – on display appears 000, the time can be adjusted for cycle 1 from 1 to 180 minutes. Pressing the button SET increases the time in minutes. After reaching the desired time, release the button SET and save it by short pressing of button START - device gives short sound signal.
7. Setting up “Cycle 2” time (Indicator Timer 2 lights) – on display appears 000, the time can be adjusted for cycle 2 from 1 to 180 minutes. Pressing the button SET increases the time in minutes. After reaching the desired time, release the button SET and save it by short pressing of button START - device gives short sound signal.

If there is no set value for Cycle 2 (000), this cycle is not executed and device works with only one cycle.



When switching on the SET UP mode temperature and time values are reset! In case of changed temperature (needed lower value) (for example from 60 to 40 degrees) it is needed to wait for the heating plate to cool down (switch off the device for at about 30 minutes).